## H130

### User Manual

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**EN**

**SAFETY NOTE**

⚠️ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.

**Introduction**

The H130 miniLab is a portable waterproof pH meter with an ISFET pH sensor for pH measurements for the food and beverage industry.

**Technical information**

|                          |   |
|--------------------------|---|
| Calibration points       | 1 point   |
| Auto buffer recognition  | pH 7.00 ; 6.86                                    |
| Temperature compensation | Automatic   |
| pH range                 | suggested range: 5 to 9 pH; range: 2 to 12 pH     |
| Temperature range        | 5 to 40 °C (40 to 105 °F)                         |
| Power                    | Auto Shut Off. 3 V lithium batteries (2x), CR2032 |

**Use the meter on**

1 Calibration

2 Measurement (direct method)

3 Cleaning

4 Battery replacement


5 Maintenance

**Precautions**


⚠️ Avoid direct contact with the pH sensor when the liquid in the reference electrode is decreased. The response time is slow or non-existent, the red status, release the electrode.

**Warranty**


1 year for manufacturing faults.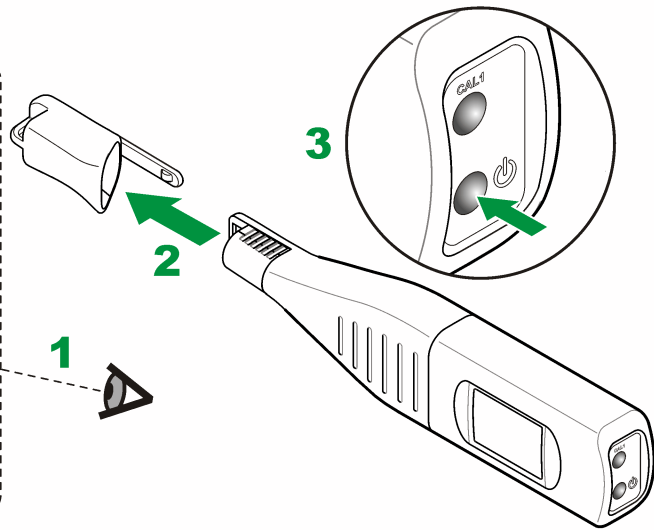


2

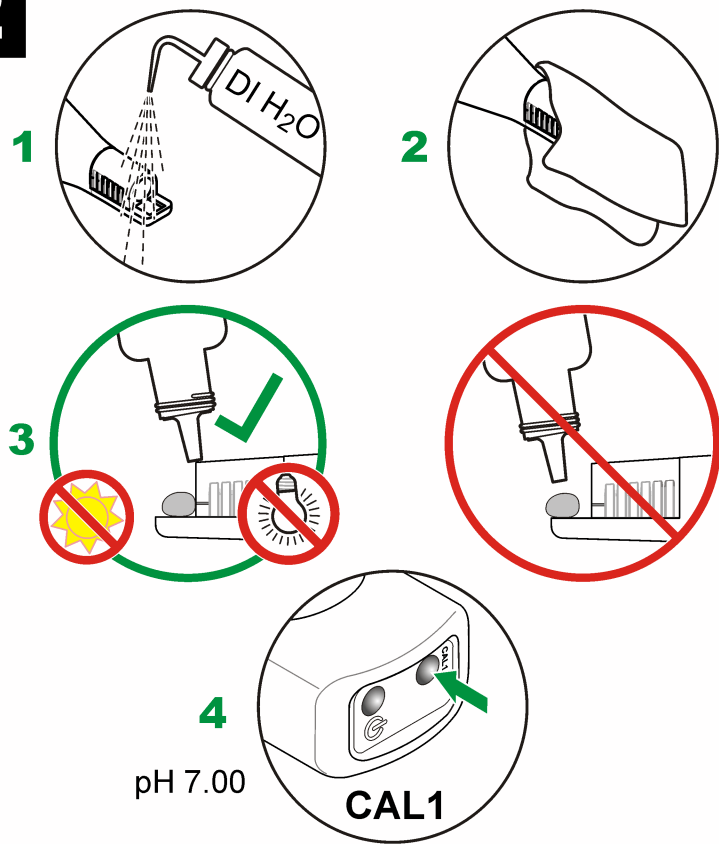
3

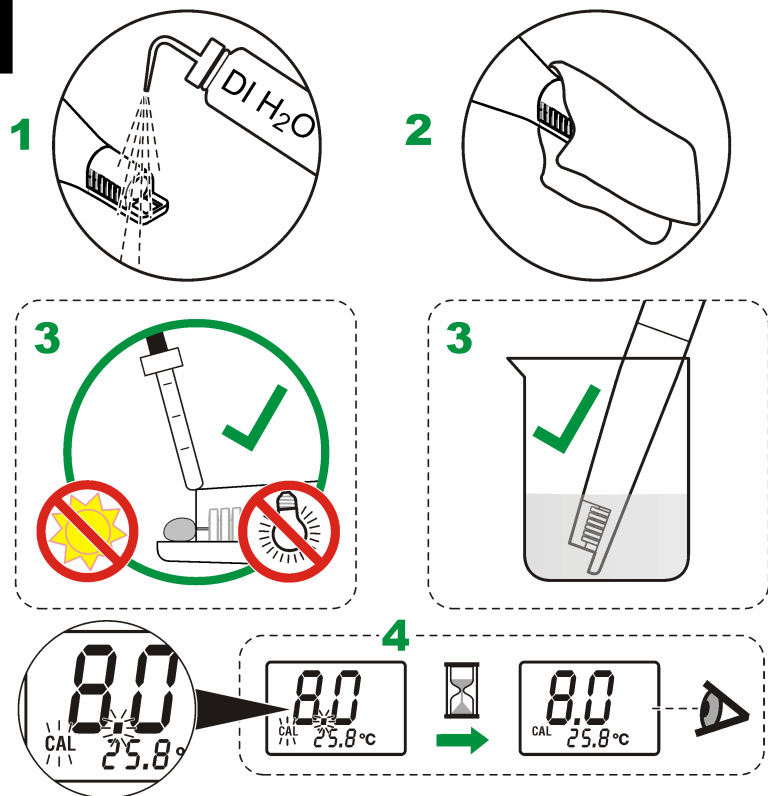
4

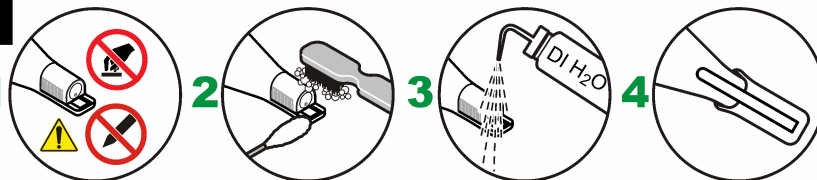5

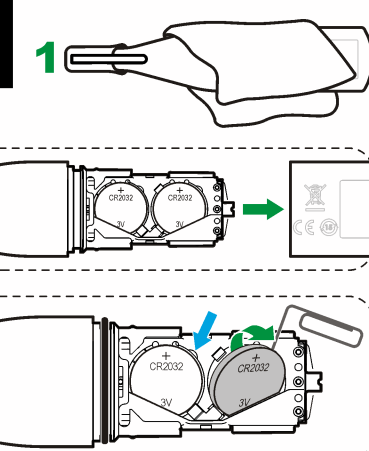6

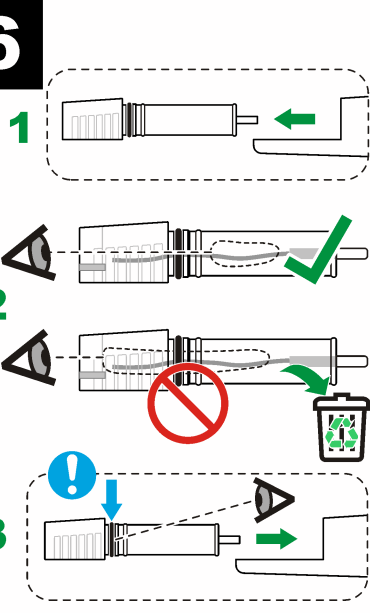EN

**Caution**

⚠️ **Chemical exposure hazard.** Refer to the current material safety data sheets (MSDS) for safety protocols.

**Introduction**

The H130 miniLab is a portable waterproof pH meter with an ISFET pH sensor for pH measurements for the food and beverage industry.

**Technical information**

|                          |   |
|--------------------------|---|
| Calibration points       | 1 point   |
| Auto buffer recognition  | pH 7.00 ; 6.86                                    |
| Temperature compensation | Automatic   |
| pH range                 | suggested range: 5 to 9 pH; range: 2 to 12 pH     |
| Temperature range        | 5 to 40 °C (40 to 105 °F)                         |
| Power                    | Auto Shut Off. 3 V lithium batteries (2x), CR2032 |

**1 Turn the meter on**

**2 Calibration**

To change the buffer set or the temperature units, push and hold CAL 1 and then push the power button. Push CAL 1 again to toggle between the parameters.

**3 Measurement**

**4 Maintenance and storage**

For protein deposits the manufacturer recommends an enzyme cleaning solution (2964349). Do not use organic solvents.

**5 Battery replacement**

Always dry the meter first.

**6 Reference electrode replacement**

Avoid direct contact with the pH sensor. When the liquid in the reference electrode is decreased, replace the electrode (9316900).

**Troubleshooting**

| Error code    | Definition   |
|---------------|--|
| Battery icon  | <10 hours battery life                             |
| BAT           | Replace the battery.                               |
| 0.00 or 14.00 | Voltage outside range. Dry or insufficient sample. |
| E06           | Slope error.                                       |
| E07           | Voltage error.                                     |
| E08           | Calibration time out.                              |
| E13           | Temperature sensor error.                          |

**Note:** When an error occurs, clean the probe and calibrate again with sufficient sample. If the problem is not solved, contact Technical Support.

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